



Daniel Kaloyanov

SpecTk 1.5.1

Facility for Rare Isotope Beams, Michigan State University, East Lansing, MI 48824 USA

2024

MICHIGAN STATE
UNIVERSITY



U.S. DEPARTMENT OF
ENERGY

Office of
Science

Changes

Grid Button

- Added a button that lets grids be applied to all selected displays

Refresh Displays

- Displays often get stuck and don't allow the user to click on displays and change them
- Added a new tool that lets the user refresh all displays on a page




Quality of Life

- Fixed the clear selected and clear page buttons
 - They no longer throw errors and will remove the data of selected displays/pages
- Added a dialog for ROI Write
 - Allows the user to choose file name and directory
- Created and implemented a icon for SpecTk
 - Allows the user to better recognize SpecTk in the taskbar

Future Updates

- Xamine Discrepancies:
 - Data index out of range
 - Counts showing differently in xamine and SpecTk

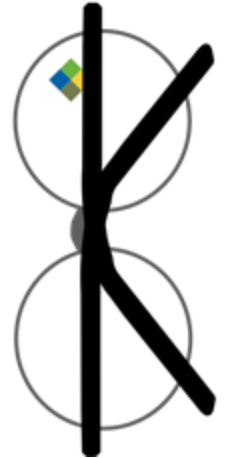

How to Use

Install/Setup

- Version 1.5.1 can be found [Here](#) or Locally at I:\projects\lisedev\Projects\SpecTclTk_Versions

Grid Button

- Select the desired displays through the following options:
 - Shift left click
 - Select Options in the bottom left of SpecTk
- Click the newly added Grid button on the left side of the screen
- The Grid button will invert the Grid value of displays even if two displays don't share Grid values

Refresh Displays

- Go to the Tool menu and select Refresh
- The tool will go line by line refreshing the displays until it runs into one of two cases
 - Finishes refreshing the page
 - Runs into a display with no Spectrum present
- To quickly fill empty displays one can use the display+ function
 - To activate it press shift while in the spectrum part of the drawer tool